



User Manual

OriCell™ Human Embryonic Stem Cell Growth Medium

Cat. No. HUXES-90011



PRODUCT DESCRIPTION:

OriCell™ Human Embryonic Stem Cell Growth Medium consists of optimized Human Embryonic Stem Cell (hESC) Basal Medium, cell culture supplements, and pre-selected fetal bovine serum. This product has been developed for the optimal expansion and maintenance of human embryonic stem cells (hESCs), thus maintaining the hESCs in the undifferentiated state while retaining the pluripotential phenotype.

The product is intended for laboratory research use only. It is not intended for diagnostic, therapeutic, clinical, household, or any other applications.

KIT COMPONENTS:

Human Embryonic Stem Cell (hESC) Basal Medium (Cat. No. HUXES-03011)	385 mL
Serum Replacement (Cat. No. SRN-19001)	60 mL
Human Embryonic Stem Cell (hESC)-Qualified Fetal Bovine Serum (Cat. No. HUXES-05001)	40 mL
Penicillin-Streptomycin	5 mL
Glutamine	5 mL
Non-essential Amino Acid	5 mL
bFGF	100 µL
2-Mercaptoethanol	500 µL

INSTRUCTIONS:

Preparation of the Complete Medium

1. Prior to use, thaw the hESC-Qualified Fetal Bovine Serum and Serum Replacement at 2-8°C overnight or until completely thawed. Gently swirl the bottles to ensure homogeneity. The serum has been heat-inactivated and is ready to use after thawing.



Note: The thawed serum may contain some flocculent precipitates. The presence of these substances in serum does not alter the performance characteristics of the product. It is not recommended to filter the serum to remove these precipitates. Doing so may result in the loss of some serum nutrients.

2. About 30 minutes prior to use, thaw Penicillin-Streptomycin solution, Glutamine solution, and Nonessential Amino Acid solution at room temperature. Gently swirl the vials to ensure homogeneity.
3. About 10 minutes prior to use, thaw bFGF and 2-Mercaptoethanol at room temperature.



Note: Centrifuge the vials briefly at low speed before removing the caps to ensure recovery of the entire content.

4. Disinfect the external surfaces of the bottles/vials for every component in the kit with 70% v/v ethanol. Allow ethanol to evaporate.
5. Aseptically open the bottles/vials inside a laminar flow hood.
6. Transfer the entire amount of hESC-Qualified Fetal Bovine Serum, Penicillin-Streptomycin solution, Glutamine solution, and Nonessential Amino Acid solution into hESC Basal Medium.
7. Rinse each vial with a small amount of basal medium. Subsequently transfer the rinse medium back into the bottle of basal medium.
8. To transfer the entire amount of bFGF and 2-Mercaptoethanol, add 0.5 mL of basal medium to the vials, mix by pipetting, and then transfer the mixture back to the bottle of basal medium as much as possible.
9. Repeat step 8 several times.
10. Gently swirl the fully supplemented (complete) medium to ensure a homogeneous mixture. The complete medium is now ready to use.



Note: Although each component in this kit is supplied sterile, it is strongly recommended to filter the fully supplemented (complete) medium.

STABILITY AND STORAGE:

All products should be stored in the dark. The hESC Basal Medium is stable at 2-8°C for up to one year. Other components are stable at -20°C for up to two years.

These products should be discarded beyond the labeled expiration date. Once prepared, the fully supplemented (complete) medium can be stored for up to one month when stored in the dark at 2-8°C.

For optimal performance, repeated warm-cooling and freeze-thawing should be avoided.

QUALITY CONTROL:

OriCell™ Human Embryonic Stem Cell Growth Medium has been tested for performance on human embryonic stem cells. The standard evaluation includes:

- Sterility test (bacteria, fungi, and mycoplasma)
- pH test
- Osmolality
- Endotoxin



Cyagen Biosciences reserves all rights on the technical documents of its OriCell™ cell culture products. No part of this document may be reproduced or adapted for other purposes without written permission from Cyagen Biosciences.